BRESSER®

**WLAN Kamera
WIFI Camera**

Art. No. 14939

DE

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EN

Instruction Manual

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About this Instruction Manual

Please read the safety instructions and the operating instructions carefully before use. Keep these instructions for renewed use at a later date. When the device is sold or given to someone else, the instruction manual must be provided to the new owner/user of the product.

General Warnings

- **Risk of choking** — Keep packaging material, like plastic bags and rubber bands, out of the reach of children, as these materials pose a choking hazard. This product contains small parts that could be swallowed by children.
- Leaking battery acid can lead to chemical burns. Avoid contact of battery acid with skin, eyes and mucous membranes. In the event of contact, rinse the affected region immediately with a plenty of water and seek medical attention.

- **Risk of electric shock** — Never bend, pinch or pull the power and connecting cables, extensions and adapters. Protect the cables from sharp edges and heat. Before operating, check the device, cables and connections for damage. Never use a damaged unit or a unit with damaged power cables. Damaged parts must be exchanged immediately by an authorised service centre.
- **Risk of property damage!** — Do not disassemble the device. In the event of a defect, please contact your dealer. The dealer will contact the Service Centre and can send the device in to be repaired, if necessary.
- Do not expose the device to high temperatures. Do not short-circuit the device or throw them into a fire. Excessive heat or improper handling could trigger a short-circuit, a fire or an explosion.
- Do not interrupt an on-going data connection to a computer before all data has been transferred. This could lead to data loss for which the manufacturer is not liable.

Introduction:

Seamlessly compatible with standard monoculars, binoculars, and astronomical telescopes, this versatile device transforms them into a powerful WIFI eyepiece camcorder. Capture stunning videos and photos from afar and store them directly in your digital camera. With 2K high-definition resolution, you can achieve clear long-distance shots. Ideal for a variety of activities such as hunting, fishing, bird watching, mountain climbing, outdoor adventures, exploration, and even concerts. Truly your ultimate companion for every adventure!

Thank you for choosing our product. Please read the user manual carefully to ensure proper installation and optimal use.

Product Description:

1. Adjustable Tightness Area
(Rotation)

2. Electronic Eyepiece

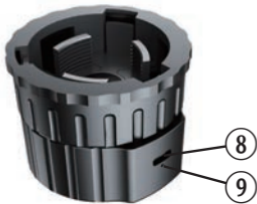
3. Fixed Buckle Position

4. TF Card Slot

5. Power On/Off Button

6. WIFI Button

7. LED Indicator light

8. USB-C Charging Port

9. Reset Hole

Installation:

Fix the camera by rotating the mounting groove to adjust the three internal buckle positions: Rotate counterclockwise to release the activity buckle. Rotate clockwise to tighten the activity buckle.

Memory Card:

Please insert the memory card before using this device. Ensure the memory card is correctly oriented with the notch prompt on the device, then press the memory card until it clicks into place. Note that the memory card must be a high-speed TF card (CLASS 10 or higher).

Power ON/OFF buttons:

Power ON: Press and hold the power button for about 3 seconds to turn on the device. The left blue light will remain on, indicating the device is on (no card status). **Start Record/Stop Record:** Short press the power button. The blue light will flash to indicate recording has started. Short press again to stop recording, and the blue light will remain on. **Power OFF:** Press and hold the power button for 3 seconds. The indicator light will turn off, indicating the device has successfully shut down.

Turn ON/OFF the WIFI:

After the eyepiece is powered on, press and hold the WIFI button for about 2 seconds. When the green light flashes, WIFI is activated and ready to connect to your mobile phone. To turn

off WIFI, press and hold the WIFI button for about 2 seconds until the green light turns off. R Reset Hole: Use this to perform a forced shutdown in case of a system crash. Simply short press the reset hole to power off the device directly.

APP Download:

Download "Onreal" App at IOS/Android APP Store in your mobile phone (or scan the QR code from the packaging box or the manual cover, as the figure shown below).

After the installation is completed, enter the APP to connect to the camera by connecting the WIFI.

WIFI name: BRESSER HD CAM

Connect WIFI and add the Device:

Here are the steps:

1. Enter the system settings WIFI on your mobile phone.
2. Locate the local WIFI hotspot named "BRESSER HD CAM" and select it to connect directly without entering a password.
3. Download the "Onreal" app from the app store.
4. Open the "Onreal" app.
5. Follow the instructions within the app to add the device.

Once the device is successfully added:

- You can configure settings and operate the camera within the app.
- You can also playback photos and videos captured by the device through the app interface.

Charge the eyepiece:

When the eyepiece is not powered on, you can charge it by inserting a USB/DC cable. The red light will turn on to indicate that charging has started. Once charging is complete, the red light will turn off. The eyepiece supports charging via USB/DC or a power bank even while it is in use.

Introduction of APP Functions:

Record: Start /Stop the recording.

Photo: Capture and save the photo

Camera File: Playback the video/photo without download, you can manage the pictures and videos from here:

Camera File

Download: Save current photo/video into Album of App, then you can view them without connect the device.

Videos screenshot: Take a picture of the currently played screen and save it as a photo.

Delete: Delete current photo/video.

Camera Settings: You can set the video resolution, sound recording on/off, speaker volume, loop duration, etc.

Camera Settings

Video Resolution: Default 2K, 2K/1080P optional.

Speaker Volume: Default Middle, Off/ Low/ Middle/ High optional.

Loop duration: 1/3/5 minutes.

Flip: Supports flipping up and down 180 degrees for video rotation.

Playback the Videos:

After downloading the video in Camera file, you can find the downloaded pictures/videos in Onreal APP Album, and then you can edit and share images and videos in the album.

File editing: Picture support image cropping, add Filter; video support video cropping and adding music.

After downloading the video, you can choose to play at a multiple speed, such as 1 frame per second or 4X multiple speed.

Multi-speed playback: When playback the video, you can choose to play at a multiple speed, such as 1 frame per second or 2X multiple speed...etc.

Overturn: Go to full screen of the video or exit full screen mode.

Connecting to PC:

After inserting a USB cable into the computer, wait a few seconds, then press the WIFI button twice. This action will automatically connect to the computer and enter USB mode.

Specifications:

CMOS/SENSOR pixel:	4MP
Video resolution:	2K@25fps/1080P@30fps
Photo resolution:	1920x1080/2304/1296
Video format:	MP4
Lens horizontal angle:	5°
Video file time:	1/3/5 minutes
WIFI point-to-point:	Support
Maximum support TF card:	128GB MAX
Battery (Li-ion)	800mAh
Recording time (WIFI) open	130 minutes
Language	Support all language photo synchronization

Notes on cleaning

- Turn off the camera
- Clean the lenses only with a soft, lint-free cloth, (e.g. microfibre cloth). To avoid scratching the lenses, use only gentle pressure with the cleaning cloth.
- Only use a dry cloth to clean the exterior of the device. To avoid damaging the electronics, do not use any cleaning fluid.

EC Declaration of Conformity

Bresser GmbH has issued a "Declaration of Conformity" in accordance with applicable guidelines and corresponding standards. The full text of the

EC declaration of conformity is available at the following internet address:

www.bresser.de/download/14939/CE/14939_CE.pdf

Disposal

Dispose of the packaging materials properly, according to their type, such as paper or cardboard. Contact your local waste-disposal service or environmental authority for information on the proper disposal.

Do not dispose of electronic devices in the household garbage!

As per the Directive 2012/19/EC of the European Parliament on waste electrical and electronic equipment and its adaptation into German law, used electronic devices must be collected separately and recycled in an environmentally friendly manner. Empty, old batteries must be disposed of at battery collection points by the consumer. You can find out more information about the disposal of devices or batteries produced after 6 January 2006 from your local waste-disposal service or environmental authority.

Drain the battery completely before disposal.

Warranty & Service

The regular warranty period is 2 years and begins on the day of purchase. To benefit from an extended voluntary warranty period as stated on the gift box, registration on our website is required.

You can consult the full warranty terms as well as information on extending the warranty period and details of our services at www.bresser.de/warranty_terms.

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Errors and technical changes reserved.

Manual_14939_WIFI-Camera_de-en_BRESSER_v072024a

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Bresser UK Ltd.